
Ref. 100150, 100152

ES.....	4	Manual original
EN.....	7	
FR.....	10	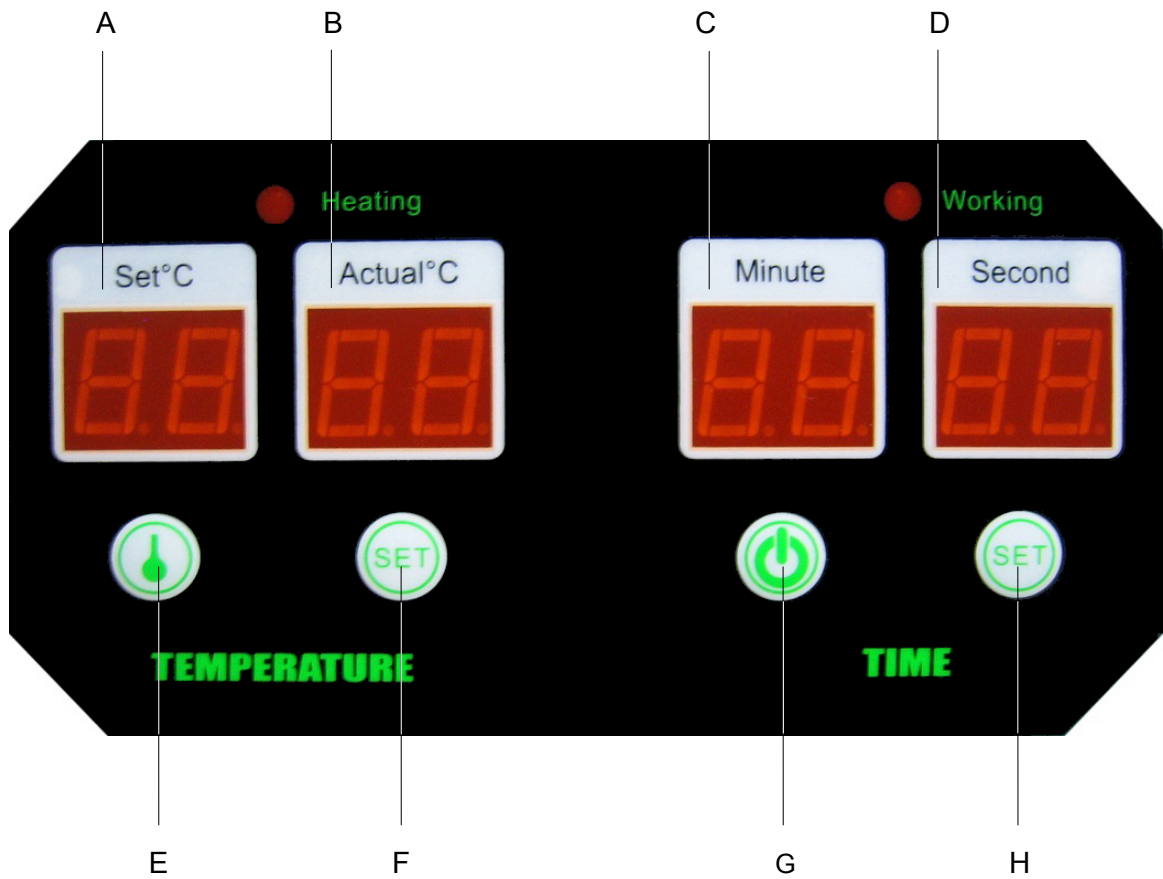

- A. Display temperatura programada / **Set temperature display** / *Display température programmée*
- B. Display temperatura real / **Actual temperature display** / *Display température réelle*
- C. Display minutos / **Minutes display** / *Display minutes*
- D. Display segundos / **Seconds display** / *Display secondes*
- E. Accionamiento temperatura / **Temperature activation** / *Activation température*
- F. Selección temperatura / **Temperature selection** / *Sélection température*
- G. Accionamiento ultrasonidos / **Ultrasound activation** / *Activation ultrasons*
- H. Selección tiempo / **Time selection** / *Sélection du temps*

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References 100150, 100152 are two practical ultrasonic baths specially designed for a wide range of applications. For a long working life and an optimum performance, we recommend you to read carefully the following instructions:

INSTALLATION

- Carefully unpack the device. Inside you will find the following accessories:
 - 1 power cord.
 - 1 plastic basket.
- Choose a flat and firm surface far from heat or vibration sources.
- Connect the machine to a grounded 230 V, 50/60 Hz power supply.
- Remove the lid, put the items in the basket. Then put the basket in the cleaning tank. In order to prevent scratches and friction of the products, do not pile them up.
- Add water or a cleaning solution in the stainless-steel tank. The liquid must cover the surface of the tools, without exceeding the fold of the tank 2 cm from the edge.
- The device is ready for operation.

OPERATION

1. Turn on the appliance main switch located at the rear.
2. Press the (H) key to set the cycle time (1 - 99 minutes):
 - Short press: increase one minute.
 - Long press: increase quickly.
3. Press the (F) key to set the operating temperature (1 - 80 °C):
 - Short press: increase one degree.
 - Long press: increase quickly.
4. Press the (G) key to start/stop the cleaning cycle.
5. Press the (E) key to activate/deactivate heating.

Do not start the appliance without liquid in the tank.

Before emptying the appliance, wait 10 minutes with the heating switched off.

MAINTENANCE AND CLEANING

To ensure optimum performance and maximum reliability of the device, we recommend:

- Do not turn on the appliance without water in the tank; The appliance may be damaged or its useful life severely shortened.
- Before emptying the appliance, wait 10 minutes with the heating switched off.
- Do not use the appliance for an excessive time or continuously.
- Do not leave liquid in the tank for a long time. After work, empty the water by opening the drain cock on the right side, clean the tank and dry it.

- Do not spill liquid on the case
- Be careful not to wet the keyboard of the device excessively.
- Use a towel to dry the tub and housing.
- Keep in a dry, cool and ventilated place.
- Do not overload the tank of the appliance with objects to be cleaned. This circumstance, in addition to reducing the effectiveness of the bath, can be detrimental to the appliance.
- Be careful not to insert excessively heavy or large objects into the tub.
- Chemical cleaning products that are compatible with the elements and parts to be treated must be used. Avoid corrosive solvents.
- When you have to clean objects for the first time - especially when they are of great value - it is advisable to carry out a test with one piece first. Ask the parts manufacturer for information on the most recommended cleaning solution.
- After using the bathroom, clean the tub and the outer casing with a damp cloth. To carry out this operation, unplug it from the network beforehand.
- To avoid short circuits, do not wash the appliance completely with water.

PRECAUTIONS

IMPORTANT- *The input voltage of the Ultrasonic Bath transducers can reach hundreds of volts. Therefore, it is very important to observe the following Safety Regulations:*

- 1) Do not open the lid when the power cord is connected to the power supply.
- 2) Do not spill water or any other conducting fluid over the unit's housing.
- 3) Always connect the unit to grounded power outlets.

- Connect the unit to a grounded 230 V, 50/60 Hz power outlet.
- Do not use inflammable solvents as cleaning solutions.
- Keep the unit unplugged before filling it. Do not fill the tank above the max. mark. Do not start up the unit when the water level is below the min. mark.
- Do not allow children or unskilled people to handle the unit.
- Never submerge the unit in any liquid to clean it.
- Ultrasonic cleaning is not suitable for organic pearls.
- Do not open the unit while it is connected to the power supply.
- The machine lid may be hot after a cycle. Proceed with caution when removing it.
- If you are not sure of the effects of the cleaning solution on a given type of item, carry out a previous test with a single item.
- Do not attempt to remove any screw or panel. The user does not need to manipulate anything inside.
- To prevent electric shock, never touch the power cord or socket with wet hands.
- Never operate the unit if the plug or the power cord is damaged. If the unit does not work properly, is damaged or has fallen into the water, refer it to a technical service for examination and repair.
- Do not touch the unit if it has fallen into the water. Unplug it immediately.
- Use the unit properly according to the user's guide. Do not use accessories not recommended by the manufacturer.
- The unit should never be left unattended when it is plugged.
- Do not operate the unit next to heat sources such as radiators.
- Install the unit on a dry and flat surface.


CLEANING SOLUTIONS

Cleaning solutions can be classified into two types: those consisting of a cleaning solution mixed with water, and solvents.


The most common cleaning method is with water, as it has an acceptable capacity to dissolve inorganic substances. However, it is extremely limited in dissolving organic substances. For this type of compound, water is a good medium for the dispersion of dirt particles, and a good basis for specific detergents that dissolve organic substances.

In most cases a mixture of water and domestic washing up liquid can be used as a cleaning solution. Nevertheless, when due to the amount or nature of dirt these cleaning solutions are not enough, specific solvents for ultrasonic baths should be used. **Under no circumstances should inflammable solvents or environmentally dangerous products be used.**


DIRT REMOVING PROCESS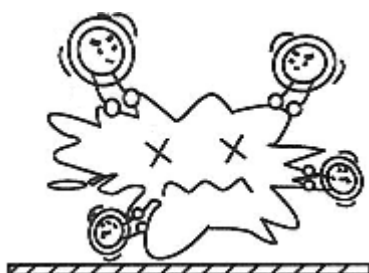

The ultrasounds generate myriads of small bubbles.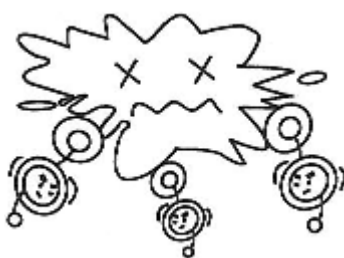

The ultrasounds make the bubbles implode constantly and impact against the object's surface

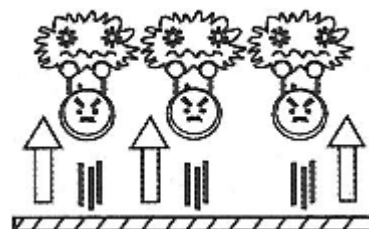The impact breaks up the dirt into many grains and separates them from the object's surface. The grains remain in the water.

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