
Ref. 080087

ES.....	9	Manual original
EN.....	10	
FR.....	11	
DE.....	12	
PT.....	13	
IT.....	14	

MESTRA[®]

DECANTADORA CON BOLSA (Tamaño reducido)
DECANTER WITH COLLECTION BAG (Small size)
BAC DE DÉCANTATION AVEC SACS (Petit modèle)
R-080087

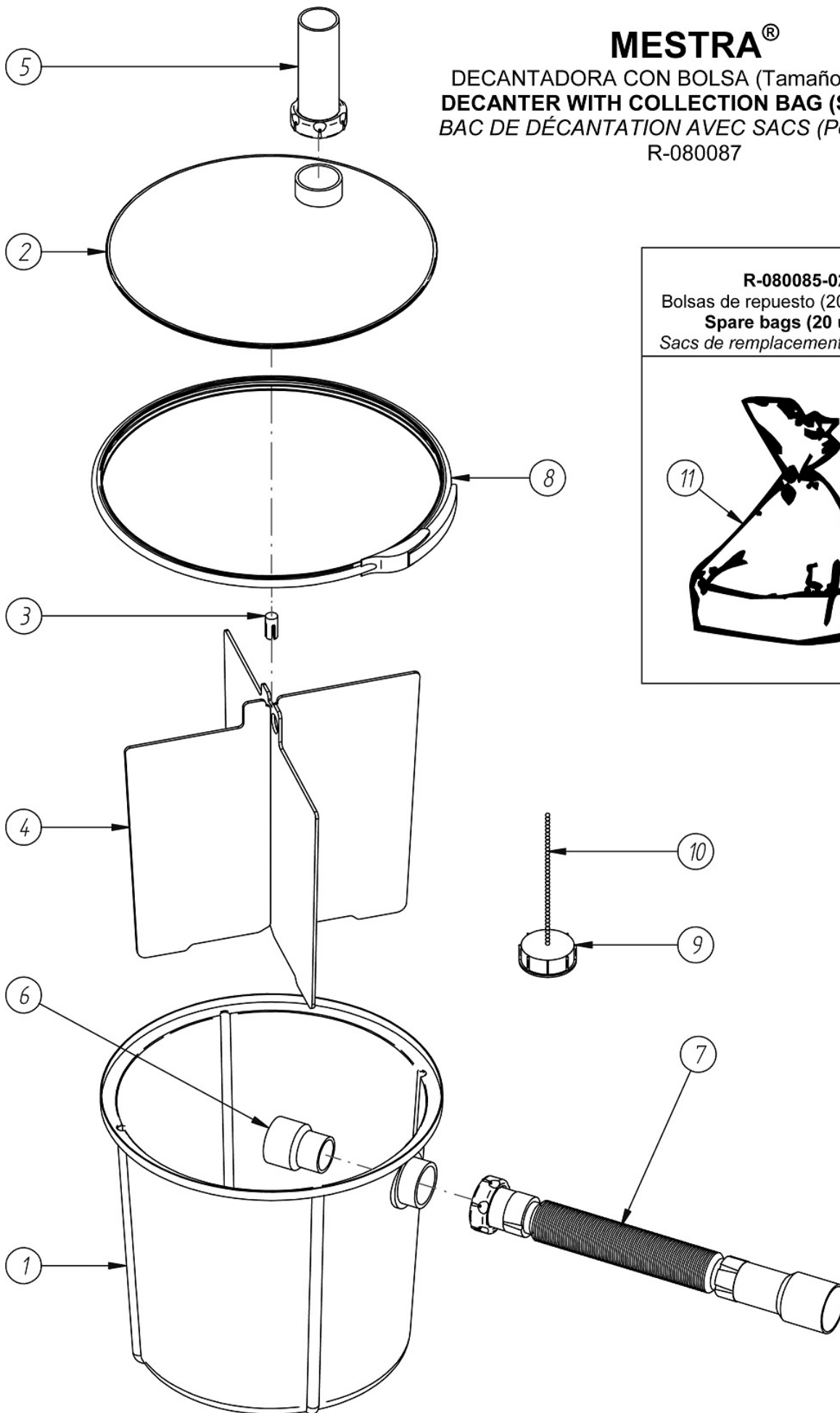
R-080085-02
Bolsas de repuesto (20 unidades)
Spare bags (20 units)
Sacs de remplacement (20 unités)

Nº	Ref.	DESCRIPTION	DESCRIPTION
1	080087-01	Cuba	Container
2	080087-02	Tapa	Lid
3	080085-08	Casquillo fijación cartabones	Squares fixation sleeve
4	080087-04	Cartabones	Squares
5	080087-05	Tubo recto entrada	Inlet straight tube
6	080087-06	Adaptador salida desagüe	Outflow exit adaptation
7	080087-07	Tubo flexible salida	Flexible exit tube
8	080087-08	Abrazadera	Hose clamp
9	080087-09	Tapón roscado	Threaded plug
10	080085-13	Cadena	Chain

Bolsas de repuesto
Spare bags
Sacs de remplacement

11	080085-02	Bolsas de repuesto (20 U.)	Spare part bags (20 U.)
----	-----------	----------------------------	-------------------------

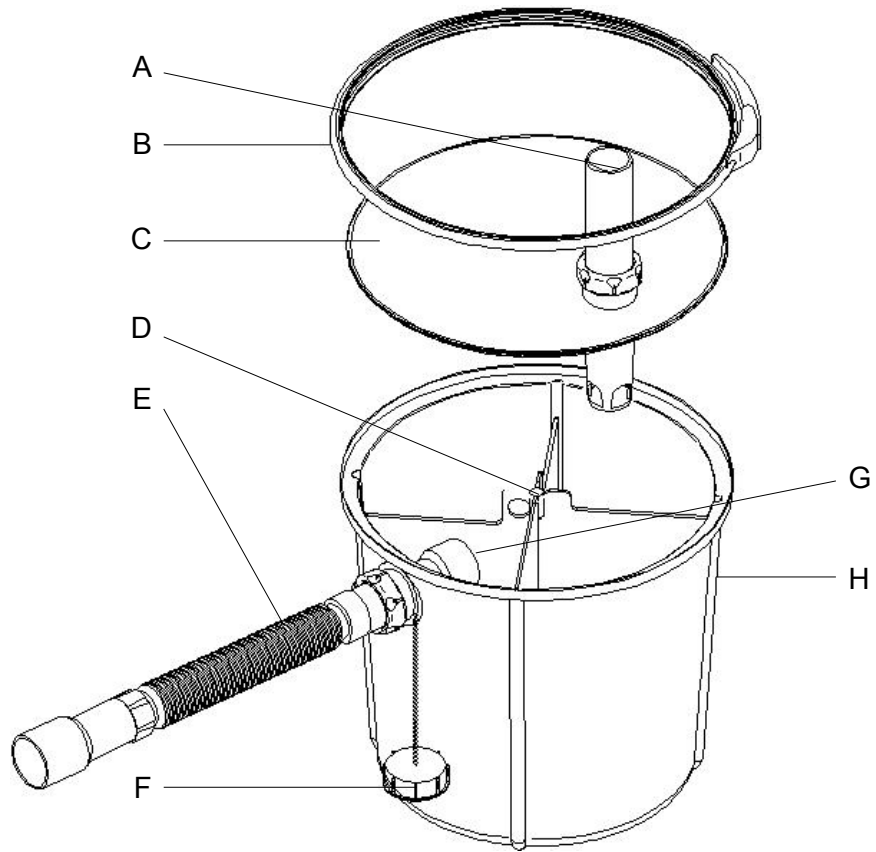ES

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CARACTERÍSTICAS TÉCNICAS

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The R-080087 decanter is made of high resistance plastic and it is specially designed to allow a quick and simple cleaning, since dirt is collected in a plastic bag. The lid closes hermetically and prevents the spread of bad odours. Its size and shape have been thought for the unit to be easily installed under the sink.

INSTALLATION (see Figures in previous pages)

1. Insert the bag into the tank while folding the bag's rim outwards.
2. Insert the separators in the grooves down to the tank bottom. The decanter's outlet should be in the lowest chamber.
3. Place the bushing on the outlet hole and press it strongly while stretching the bag.
4. Open the drain hole in the bag with a knife.
5. Once the drain hole has been opened, press the bushing until it remains correctly fitted.
6. Place the lid making sure the water inlet is in the highest chamber.

In order to ensure an optimum operation of the decanter and to prevent deformations, the lid should be closed observing the following instructions:

7. Place the decanter's lid over the tank and make sure it is well centred and fitted all along its perimeter. Open the clamp completely and place it around the decanter.
8. Loosen the clamp's clasp and make sure the lid is well centred and fitted before tightening the clamp.
9. Tighten the clamp carefully to prevent the lid moving.
10. Close the clamp completely. The lid should be perfectly fitted to the tank.
11. Put the decanter under the sink. Connect the sink's outlet to the decanter's inlet and the latter's outlet to a drainpipe. In case of any trouble when adapting the decanter to the drainage system in your lab, consult a professional plumber.

CLEANING

12. To remove the decanter from under the sink, loosen the inlet and outlet threads.
13. Put the threaded plug on to close the drain hole. Open the lid and remove the remaining water.
14. Remove the separators.
15. Remove the adaptor from the outlet hole.
16. Tie up the bag below the drain hole and dispose of the waste.

NOTE: when cleaning the decanter, check that there are no obstructions in the drain pipes.

CAUTION

- Place the decanter on a firm, flat and balanced surface.
- Make sure there are not electric wires or circuits near the decanter. A minimum safety distance of 1 meter is recommended.
- Be extremely cautious while transporting or moving the decanter. Prevent the risk of turning the unit over.
- In case of any trouble when adapting the decanter to the drainage system in your lab, consult a professional plumber.

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