
R-050051/52/53/54

ES..... MECHERO DE GAS	3
EN GAS BURNER.....	5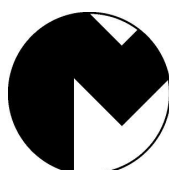

MESTRA®


TALLERES MESTRAITUA S.L.

Txori-Erri Etorbidea, 60

Tel. (+34) 944530388 - Fax (+34) 944711725

mestra@mestra.es - www.mestra.es

48150 SONDIKA - BILBAO - ESPAÑA

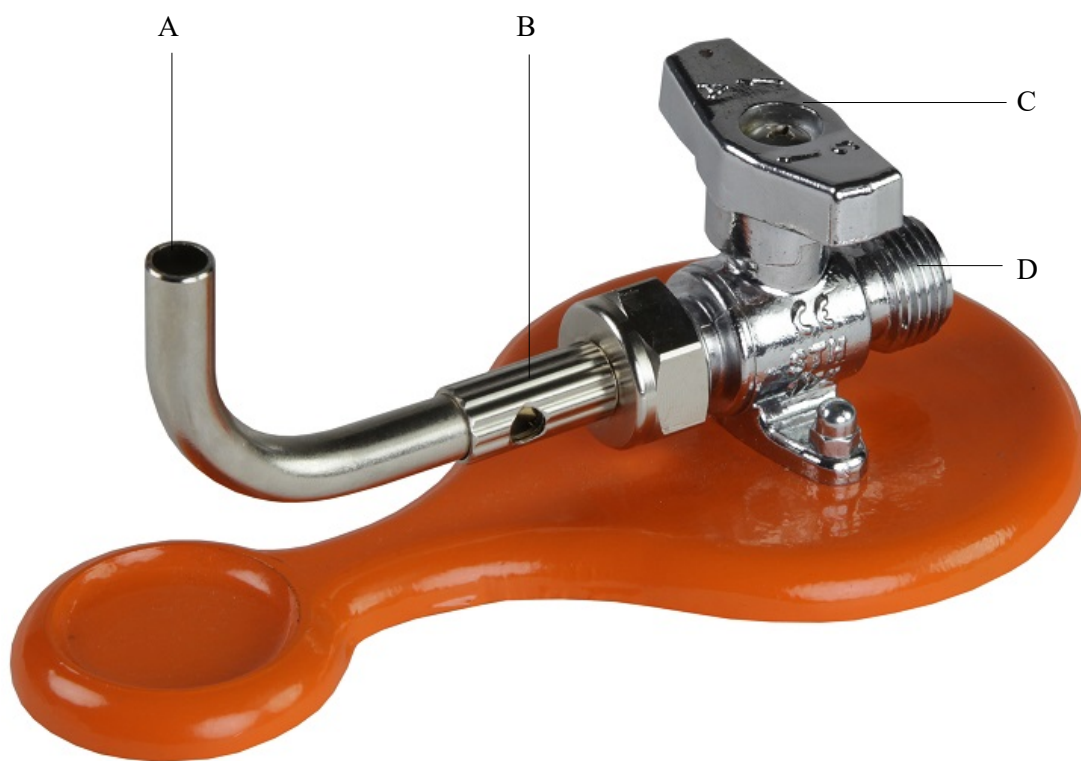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EN

INSTALLATION

- Carefully unpack the gas burner. You will find a label on the rear side of the burner base describing the type of fuel to be used. **Make sure that the burner fuel coincides with your gas supply.**

R-050051 Butane/propane gas

R-050052 Natural gas

R-050053 Butane/propane gas + wax container

R-050054 Natural gas + wax container

- Place the burner on a flat surface with no vibrations at a well-ventilated area, away from flammable materials.
- Check your gas supply conditions. Make sure that all the necessary inspections have been performed by licensed technicians.
- To connect the burner to the gas supply, screw a normalized hose to the burner fitting (D). Make sure that the nut is correctly tightened.

In case you are using a connection dowel (look at the "Accessories" section), insert the dowel into the rubber tube and tighten it with a clamp. Then, screw the dosel to the burner fitting (D).

USE OF THE BURNER

1. If you need, place the wax container (look at the “Accessories” section) upon the flame nozzle (A).
2. Turn the air flow regulator (B) until the circular hole is completely open.

3. Open the main stopcock of your gas supply.
4. Then, press the knob of the gas tap (C) downwards in order to unlock. Then apply a ¼ turn until it is parallel to the central axis of the burner.

Close

Open

5. Bring up a flame to the nozzle (A). The burner should immediately light up with no problem.
6. You can use the regulator (B) to adjust the airflow. By opening the regulator completely, you will obtain a flame with a blue base and a bright yellow tip. Length and bright will be constant with no alterations. This kind of flame means that combustion is perfect and the gas supply is optimum.

If you use the regulator (B) to shut the airflow off, the blue color will eventually become reddish, meaning that combustion is lacking oxygen.

Under extreme circumstances, when the airflow is too low, black smoke (unburned) may appear. **This situation should always be avoided, since the resulting gases are extremely dangerous.**

7. When work is finished, close the gas tap (C) applying a ¼ turn until it is perpendicular to the burner axis. Then close the gas main stopcock.

MAINTENANCE AND CLEANING

1. Gas facilities must be inspected regularly by licensed technicians.
2. Clean the burner externally on a regular basis. First make sure that the gas mains stopcock has been shut off completely. Use a cloth soaked in soapy water to clean the unit. Never use alcohol or flammable products.
3. During normal operation, molten wax may accidentally fall through the orifice in the flame nozzle (A). Wax accumulating inside the nozzle elbow may cause obstructions and impair burner performance. Should this occur, proceed to remove the wax as follows:
 - Close the main stopcock.
 - Release the nut of the hose from the connection fitting (D).
 - **Once the burner has been disengaged from the main supply**, apply a ¼ anti-clockwise turn to the flame nozzle (A).
 - Apply a flame to the nozzle elbow to melt the wax accumulated inside. Tilt the burner to allow the wax to flow out freely through the flame nozzle (A).

- After removing the wax, apply a ¼ turn to the flame nozzle to return to the initial position, and connect the burner to the gas mains as described above, in reverse order.

PRECAUTIONS

- Before connecting the burner, ensure that the gas supply is correct (butane, natural gas).
- Only connect the appliance to a gas network that has been subjected to regulatory inspections by an accredited technician.
- Make sure that the connection between burner and gas supply is solid and secure.
- Never allow children or unqualified staff to use the burner.
- Make sure the flame has a bright blue base and a strong yellow tip. Reddish colors, and black smoke in particular, indicate poor combustion. Always avoid this situation using the airflow valve.
- After using the burner, always close the shutoff cock and gas mains stopcock.
- Place the burner on a flat surface with no vibrations at a well ventilated area, away from flammable materials.
- Never leave the burner operating unattended.
- If you smell gas, close the mains stopcock immediately and inform a licensed technician.
- Make sure the supply hose is always in perfect condition as required by regulations, as it could cause an explosion or Accidental poisonings. Check the expiration date and, if the rubber is expired, is deformed or worn, it is imperative to switch to another of the same type and length in good condition.

ACCESSORIES

- 050050-01 Wax container:
Burner series R-050053/54 is supplied with a wax container.
Burner series R-050051/52, is identical to the previous one, but without this accessory.
In order to use the container, place it upon the flame nozzle.

- 050051-04 Connection dowel:
Components:

Assembly:

TECHNICAL CHARACTERISTICS

	050051/050053	050052/050054
Gas:	G30 (Butane) / G31 (Propane)	G20 (Natural)
Category:	I _{2H}	I ₃₊
Pressure:	28-30 / 37 mbar	20 mbar
Injector:	0.3 mm	0.5 mm
Consumption:	Hi: 0.3 kW	
Height:	72 mm	
Length:	175 mm	
Width:	94 mm	