
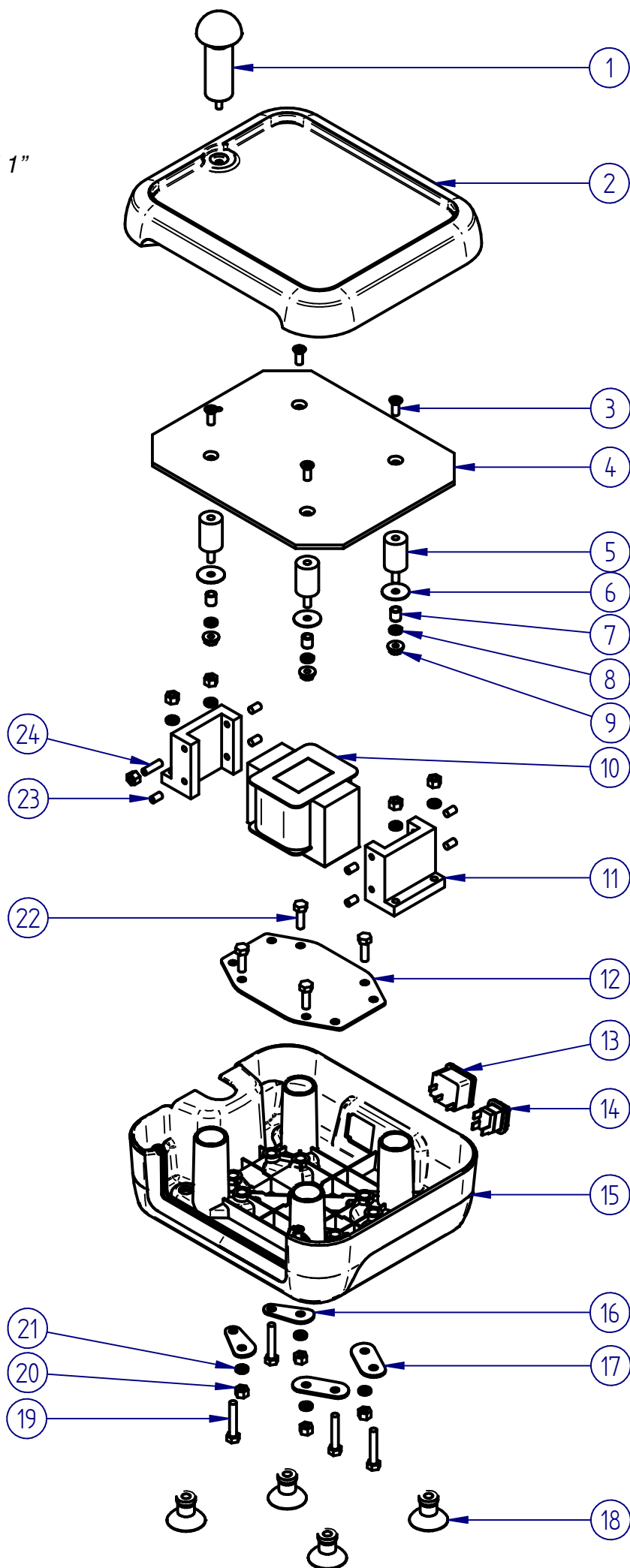
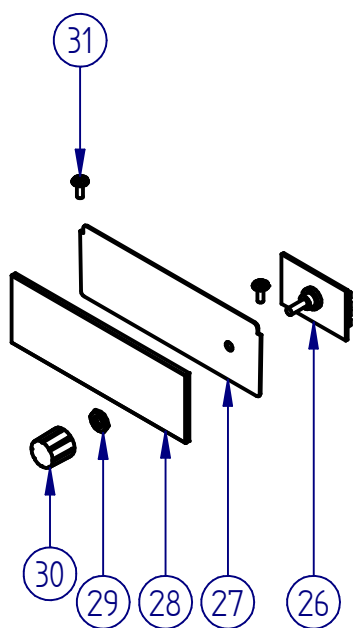


# MESTRA<sup>®</sup>

ELECTRO VIBRADOR "TEIDE 1"  
ELECTROVIBRATOR "TEIDE 1"  
VIBRATEUR ELECTRIQUE "TEIDE 1"  
R-080021

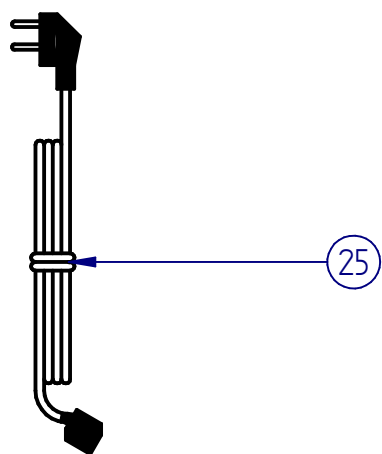

# MESTRA®

ELECTRO VIBRADOR "TEIDE 1"  
**ELECTROVIBRATOR "TEIDE 1"**  
VIBRATEUR ELECTRIQUE "TEIDE 1"  
Ref. 080021

Nº	Ref.	DESCRIPCIÓN	DESCRIPTION
1	080022-10	Bola de vaciado	Emptying ball
2	080022-06	Plataforma de goma	Rubber platform
3	080020-34	Tornillo plataforma	Platform bolt
4	080022-07	Plataforma metálica	Metal platform
5	080020-31	Silent block	Silent block
6	080022-17	Arandela grande silent block	Silent block large washer
7	080021-03	Casquillo	Bushing
8	040064-62	Arandela pequeña silent block	Silent block small washer
9	080022-18	Tuerca con arandela	Nut with washer
10	080020-11	Electroimán 230 V	Electromagnet 230 V
	080020-11B	Electroimán 115 V	Electromagnet 115 V
11	080020-14	Soporte electroimán	Electromagnet support
12	080022-20	Soporte base	Base support
13	080490-06	Base de conexión	Connection base
14	080422-17	Interruptor general	Main switch
15	080022-03	Carcasa	Housing
16	080022-22	Soporte exterior pequeño	Small out support
17	080022-21	Soporte exterior grande	Large out support
18	080020-30	Ventosa	Suction cap
19	040064-32	Tornillo electroimán	Electromagnet bolt
20	080022-15	Tuerca	Nut
21	080022-14	Arandela	Washer
22	060064-28	Tornillo soporte	Support bolt
23	080020-15	Espárrago corto	Short stud
24	080010-16	Espárrago largo	Long stud
25	000250	Cable 230 V	Cable 230 V
	000252	Cable 115 V	Cable 115 V
26	080030-12	Tarjeta control	Control board
27	080021-02	Soporte control	Control support
28	080021-01	Carátula control	Control cover
29	080010-18	Tuerca control	Control nut
30	080010-19	Pomo control	Control knob
31	080093-32	Tornillo control	Control bolt